



ECOPLUS

60X-100X MICROSCOPE

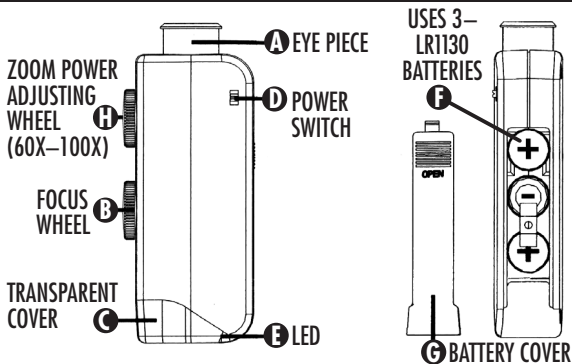


TO OPERATE MICROSCOPE

- 1) Place specimen on flat surface.
- 2) Place microscope upright with transparent cover (C) face down.
- 3) Switch on LED (E) by sliding Power Switch (D).
- 4) Adjust Zoom Power Adjusting Wheel (H) for magnifying power (60X-100X).
- 5) Look through eyepiece (A), turn Focusing Wheel (B) to obtain sharpest image.

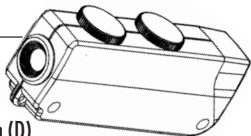

TO INSTALL BATTERIES

- 1) Slide out battery cover (G) from rear.
- 2) Install 3-LR1130 (F) batteries with (+) positive at either end & (-) negative in the middle.
- 3) Slide Battery Cover (G) firmly back into place & the microscope is ready to use.

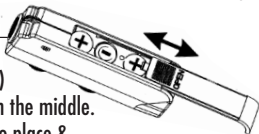

NOTE: Replacement of LED not recommended